



Lifespan Technologies
PO Box 711297
Salt Lake City, UT 84171
Tel: 801-971-2020 Fax: 801-464-6116
info@lifespantech.com

Product No: **K-5800**

K-5800 Dermatan sulfate / Heparan
Sulfate ELISA for buffer

K-5800: Dermatan / Heparan Sulfate ELISA for Buffer or Urine Samples

INTENDED USE: THIS PRODUCT IS FOR RESEARCH USE ONLY. NOT INTENDED FOR CLINICAL OR DIAGNOSTIC USE.

Kit includes:

Coated 96-well plate
Detector -Enzyme Conjugate vial
Conjugate Diluent
TMB Solution
Stop Solution, 0.5M H₂SO₄
Wash Concentrate 10X, (dilute 1 part plus 9 parts water to make TBS plus 0.05% TWEEN 20)


Researcher must provide:

Pipettes (8-channel multipipettor recommended)
Absorbance microplate reader
Dermatan / heparan sulfate standards
Normal Saline (154 mM NaCl)
Plate Cover

Storage and Stability

Kit can be stored unopened at 4°C for up to six months. Opened and reconstituted solutions can be used for up to one week when stored at 4°C. All components and solutions should be protected from light. Reconstituted detector enzyme conjugate should be used immediately or aliquotted and stored at -80°C.

Background

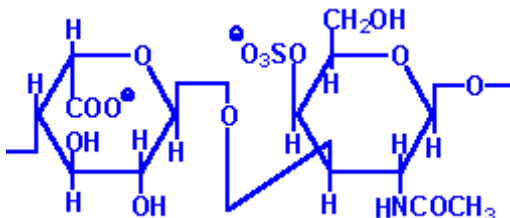

Dermatan sulfate (DS) is in the family glycosaminoglycan. DS is found mostly in the skin but also in blood vessels, tendons, heart valves, and pulmonary connective tissue. Heparan sulfate (HS) is a sulfated polysaccharide that is found on the surface of most cells as part of proteoglycans. Heparan sulfate is also present in the extracellular matrix. The polysaccharide mediates the interactions between a number of different proteins.

The DS / HS ELISA, product number K-5800, is a quantitative enzyme-linked assay designed for the *in vitro* measurement DS / HS levels in buffer. This assay measures DS / HS directly using a DS / HS binding protein which has been conjugated to HRP.


The DS / HS -ELISA is a competitive assay in which the colorimetric signal is inversely proportional to the amount of DS / HS present in the sample. Samples to be assayed are first mixed with the detector-enzyme conjugate in wells of the coated plate. DS / HS in the sample competes with DS / HS bound to the plate for binding of the detector-enzyme conjugate. The concentration of DS / HS in the sample is determined using a standard curve of known amounts of DS / HS.

Reagent Preparation

Dermatan / Heparan Sulfate Standards: Prepare standards using fresh normal human plasma and your DS / HS to obtain standards of 0.03, 0.1, 0.3, 1.0, 3.0 and 10.0 µg/mL. **Standardization should be performed using DS / HS that is the same DS / HS type contained in your unknowns.**

Working Detector-Enzyme Conjugate: Measure exactly **9.1mL** of conjugate diluent and add to a clean tube. Perform a 'clean transfer' of the tube of the Detector-Enzyme Conjugate into the 9.1 mL of conjugate diluent. This can be done by adding 500 microliters of the measured Diluent to reconstitute the Detector-Enzyme. Wait a minute to allow the lyophilized material to dissolve and then add the liquid back to the tube. Repeat this step two more times to be sure all the Detector-Enzyme Conjugate has been transferred from the vial to the tube. Reconstituted Detector-Enzyme Conjugate is unstable in the Conjugate Diluent and should be used immediately. If you wish to perform less than a full plate, the Reconstituted Detector-Enzyme Conjugate can be stored frozen at -80°C and thawed immediately before re-use.

1X Wash Buffer: Make a 1:10 dilution of 10X Wash B₁

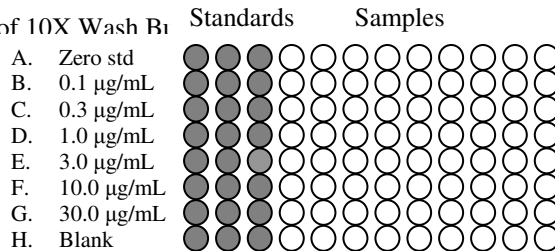

Dermatan/Heparan Sulfate Plate

Assay Procedure

1. Set up the ELISA plate as illustrated above. We suggest the standard dilution series be run in triplicate for best results. Add **50 µL** of Standards and samples into corresponding wells. Add **50 µL** of Working Detector-Enzyme Conjugate to all wells except the Blank wells. Mix well. Cover plate and incubate for one hour at room temperature. A rotator is recommended if available.
2. Discard the solution and wash the wells four times with 300 µL per well of 1X Wash Buffer. An automated plate washer is recommended if available. After washing, immediately proceed to the next step. Do not delay in removing wash buffer from the wells. Do not allow plate to dry.
3. Add 100 µL TMB Solution to each well. Incubate the plate in the dark at room temperature for 40-60 minutes waiting for the zero wells to develop to a medium to dark blue color. Watch for color development and DO NOT overdevelop.
4. Add 50 µL Stop solution which will change the color from blue to yellow.
5. Immediately measure the absorbance of each well at 450 nm.
6. Calculate the binding percentage for each sample using the formula:

$$[A_{450}(\text{Sample}) - A_{450}(\text{Blank})] / [A_{450}(\text{Zero Dermatan Sulfate}) - A_{450}(\text{Blank})] \times 100 = \% \text{ Binding}$$

Using linear or nonlinear regression, plot a standard curve of percent binding versus concentration of standards. Determine levels of unknowns by comparing their percentage of binding relative to the standard curve. Concentrations can be estimated by comparing the values from the wells containing unknowns to the values in the standard curve.

Lifespan Technologies products are sold for research and development purposes only and are not for diagnostic use or to be incorporated into products for resale without written permission from Lifespan Technologies. Materials in this publication, as well as applications, methods and use, may be covered by one or more U.S. or foreign patents or patents pending. We welcome inquiries about licensing the use of our trademarks and technologies at info@lifespantech.com.