



Lifespan Technologies
PO Box 3352
Salt Lake City, UT 84110
Tel: 801-971-1818 Fax: 801-464-6116
info@lifespantech.com

Product No: **K-2300**

Unfractionated Heparin (UFH)
Ultra Range (0.3 U/ml – 100 U/mL)

Unfractionated Heparin (UFH) Ultra Range ELISA Kit for Buffer Samples

INTENDED USE: THIS PRODUCT IS FOR RESEARCH USE ONLY. NOT INTENDED FOR CLINICAL OR DIAGNOSTIC USE.

Kit includes:

Heparin-coated 96-well plate
Detector -Enzyme Conjugate
Conjugate Diluent
TMB Solution
Stop Solution
Wash Concentrate 10X, (dilute 1 part plus 9 parts dH₂O)


Researcher must provide:

Multichannel Pipetter
Absorbance microplate reader
Plate Cover
UFH Standard from USP reference or your heparin
Standard Diluent, Tris Buffered Saline (TBS) pH 7.5 (10mM Tris, 150mM NaCl)

Storage and Stability

Kit can be stored unopened at 4°C for up to six months. Reconstituted detector enzyme conjugate is unstable and should be used immediately. If you wish to run less than a full plate, it should be stored as frozen aliquots at -80°C. Aliquots must be thawed **immediately** before use. After one thaw, any unused detector enzyme conjugate must be discarded. The TMB solution should be protected from light.

Background

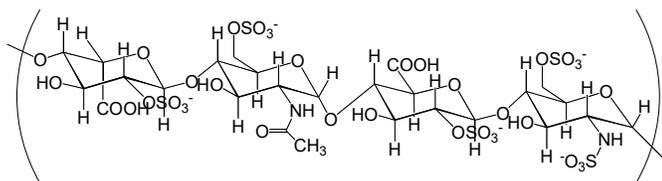

Heparin (heterogeneous)

Heparin is a glycosaminoglycan with alternating uronic acid and aminoglycoside units. It is an anticoagulant used either in its native unfractionated form (UFH) MW ~16 kD or in various partially depolymerized forms (LMWH) of 4-8 kD. The heparin-ELISA product number K-2300 is a quantitative enzyme-linked assay designed for the *in vitro* measurement of unfractionated heparin levels in low protein content fluids such as buffer or urine. This

assay measures heparin directly using a heparin binding protein which has been conjugated to HRP.

The heparin-ELISA is a competitive assay in which the colorimetric signal is inversely proportional to the amount of heparin present in the sample. Samples to be assayed are first mixed with the Detector-Enzyme Conjugate in wells of the heparin coated plate. Heparin in the sample competes with heparin bound to the plate for binding of the Detector-Enzyme Conjugate. The concentration of heparin in the sample is determined using a standard curve of known amounts of heparin. **For best results, the heparin used for the standard curve should match the type of heparin being assayed.**


Reagent Preparation

Heparin Standards: Make dilutions of your Heparin Standard using Tris Buffered Saline (TBS) pH 7.5 (10mM Tris, 150mM NaCl) to obtain standards of 0.3, 1.0, 3.0, 10.0, 30, 100 U/mL. **Standardization should be performed using heparin that is the same heparin type contained in your unknowns.**

Working Detector-Enzyme Conjugate: Measure exactly **9.1 mL** of conjugate diluent and add to a clean tube. Perform a 'clean transfer' of the lyophilized Detector-Enzyme Conjugate into the 9.1 mL of conjugate diluent. This can be done by adding 500 microliters of the measured Diluent to reconstitute the Detector-Enzyme. Wait a minute to allow the lyophilized material to dissolve

and then add the liquid back to the tube containing the conjugate diluent. Repeat this step two more times to be sure all the Detector-Enzyme Conjugate has been transferred from the vial to the tube. If you wish to perform less than a full plate, the reconstituted Detector-Enzyme Conjugate must be stored as aliquots at -80°C. Aliquots must be thawed immediately before use. After one thaw, any unused detector enzyme conjugate must be discarded.

1X Wash Buffer: Make a 1:10 dilution of 10X Wash Buffer in distilled or deionized water.

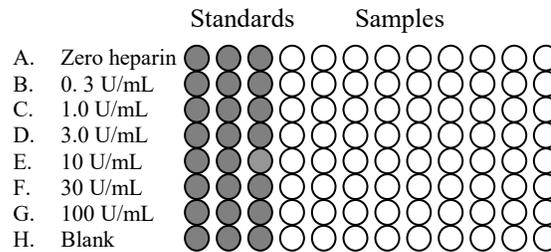

Heparin ELISA Plate

Assay Procedure

1. Set up the heparin ELISA plate as illustrated above. We suggest the heparin standard dilution series be run in triplicate for best results. Add **10 µL** of Standards and samples into corresponding wells. Add **90 µL** of Working Detector -Enzyme Conjugate to all wells except the Blank wells. Mix well by gently tapping plate from side to side. Cover plate and incubate for 30 minutes at room temperature. A rotator is highly recommended if available, as constant mixing significantly improves precision.
2. Discard the solution and wash the wells four times with 300 µL per well of 1X Wash Buffer. An automated plate washer is recommended if available to improve precision. After washing, immediately proceed to the next step. Do not delay in removing wash buffer from the wells. Do not allow plate to dry.
3. Add 100 µL TMB Solution to each well. To improve precision, do not pipette bubbles into wells. Incubate the plate in the dark at room temperature for 4-60 minutes waiting for the zero heparin wells to develop to a medium to dark blue color. Periodically check for color development and **DO NOT** overdevelop.
4. Add 50 µL Stop Solution which will change the color from blue to yellow.
5. Immediately measure the absorbance of each well at 450 nm using the absorbance microplate reader.
6. Calculate the binding percentage for each sample using the formula:

$$[A_{450}(\text{Sample}) - A_{450}(\text{Blank})] / [A_{450}(\text{Zero heparin}) - A_{450}(\text{Blank})] \times 100 = \% \text{ Binding}$$

Using linear or nonlinear regression, plot a standard curve of percent binding versus concentration of heparin standards. Determine heparin levels of unknowns by comparing their percentage of binding relative to the standard curve. Heparin can be estimated by comparing the values from the wells containing unknowns to the values in the standard curve.

Lifespan Technologies products are sold for research and development purposes only and are not for diagnostic use or to be incorporated into products for resale without written permission from Lifespan Technologies. Materials in this publication, as well as applications and methods and use, may be covered by one or more U.S. or foreign patents or patents pending. We welcome inquiries about licensing the use of our trademarks and technologies at MM@lifespantech.com.